DOS G 80 flex  
DOS G 80/1 flex  
DOS G 80 ProfiLine flex  
DOS G 80/72

<b>en</b>	Fitting instructions Dispensing module
<b>it</b>	Montavimo instrukcija DOS modulis
<b>lv</b>	Montāžas instrukcija DOS modulis
<b>nl</b>	Montagehandleiding DOS-module
<b>no</b>	Monteringsveiledning DOS-modul
<b>pl</b>	Instrukcja montażu Moduł DOS
<b>pt</b>	Instruções de montagem Módulo DOS
<b>ru</b>	Инструкция по монтажу DOS-модуль
<b>sk</b>	Návod na montáž DOS modul
<b>sl</b>	Navodila za montažo Dozirni modul
<b>sv</b>	Monteringsanvisning DOS-modul
<b>tr</b>	Montaj Talimatı DOS Modülü
<b>zh</b>	安装说明 配给模块

**en** ..... 4  
**lt** ..... 14  
**lv** ..... 24  
**nl** ..... 34  
**no** ..... 44  
**pl** ..... 54  
**pt** ..... 64  
**ru** ..... 74  
**sk** ..... 87  
**sl** ..... 97  
**sv** .....107  
**tr** .....117  
**zh** .....127

---

<b>Product description</b> .....	4
<b>Caring for the environment</b> .....	6
Disposal of the packing material .....	6
Disposing of your old machine.....	6
<b>Installation</b> .....	7
Schematic diagram showing the dispensing module .....	7
Fitting the dispensing module .....	7
Wall mounted: .....	7
Surface mounted:.....	7
Installation inside an adjacent unit: .....	8
<b>Connection</b> .....	9
Connecting the dispensing modules.....	9
DOS G 80 with hose adapter .....	9
DOS G 80 ProfiLine with hose adapter and dispenser connection .....	9
Strain relief for dispensing hoses .....	10
Connecting the dispensing hoses.....	10
Maintenance.....	11
<b>Technical data</b> .....	12

## en - Product description

---

The dispensing module is an external dispensing system that is intended solely for connection to Miele cleaning machines. It dispenses liquid process chemicals (detergent or rinsing agent, for example) automatically at the appropriate stage of a programme. The dispensing module does away with the need to fill the dispenser containers in the door of the machine (if there are any). Up to two external dispensing systems can be connected depending on the type of machine.

Suitable process chemicals are available from Miele Professional.

Dispensing modules can be retrofitted and connected if required.

 Risk posed by conversions or modifications

Please note that the dispensing module must only be used for the applications described in these instructions.

Any other use, conversion or modification is not permitted and potentially dangerous.

Miele cannot be held liable for damage resulting from incorrect or improper use or operation.

Initial commissioning will be undertaken by the Miele Customer Service Department or authorised service technicians, who will provide you with operating instructions.

Please follow the operating instructions for your cleaning machine and pay particular attention to warning and safety instructions when using and handling chemical agents.

 Damage caused by noncompliance with the fitting instructions.

Please read these fitting instructions carefully before fitting the dispensing module.

This prevents both personal injury and damage to the dispensing module.

Keep these fitting instructions in a safe place.

### Areas of application for dispensing modules

Dispensing modules are suitable for use with the following commercial dishwashers and washer-disinfectors:

#### DOS G 80

- |              |             |           |
|--------------|-------------|-----------|
| – PG 8055    | – PG 8060   | – PG 8096 |
| – PG 8056    | – PG 8061   | – PG 8099 |
| – PG 8057 TD | – PG 8063   | – PG 8504 |
| – PG 8058    | – PG 8067 * | – PG 8583 |
| – PG 8059    | – G 8066 *  | – PG 8593 |

\* When connecting the dispensing module to these commercial dishwashers please use the hose adapter supplied and double the dispensing concentration.

#### DOS G 80 ProfiLine

- |                  |                  |
|------------------|------------------|
| – PG 8080 *      | – PG 8130 *      |
| – PG 8081i *     | – PG 8131i *     |
| – PG 8082 SCi *  | – PG 8132 SCi *  |
| – PG 8083 SCVi * | – PG 8133 SCVi * |

\* When connecting the dispensing module to these commercial dishwashers please use the hose adapter supplied and fit the dispenser connection.  
For PG 8080, PG 8081i, PG 8082 SCi and PG 8083 SCVi commercial dishwashers, it is also necessary to double the dispensing concentration.

#### DOS G 80/72

- PG 8172

### Symbols on the DOS module

Liquid agent flow direction

Warning:  
Danger of electric shock


### **Disposal of the packing material**

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in landfill sites.

### **Disposing of your old machine**


Electrical and electronic machines often contain valuable materials. They also contain materials which, if handled or disposed of incorrectly, could be potentially hazardous to human health and to the environment. They are, however, essential for the correct functioning of your machine. Please do not therefore dispose of it with your household waste.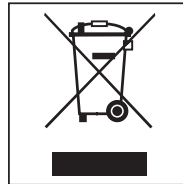

Please dispose of it at your local community waste collection / recycling centre or contact your dealer for advice.

Ensure that it presents no danger to children while being stored for disposal.

## Schematic diagram showing the dispensing module

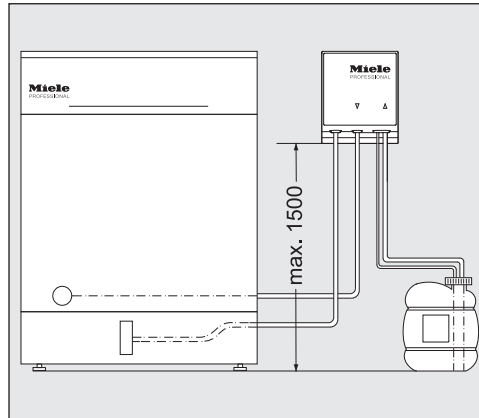
⚠ Dispensing problems caused by fitting the canister incorrectly. The canister must not be placed on top of or above the machine. Make sure that the dispensing hose is not kinked or trapped. Place the canister on the floor next to the machine or inside an adjacent unit.

## Fitting the dispensing module

The dispensing module is supplied with two screws (4.2 x 25 mm) and two plugs (6 mm) for fitting.

- Select a suitably accessible position, e.g. on an adjacent wall or inside a unit, for fitting the dispensing module.

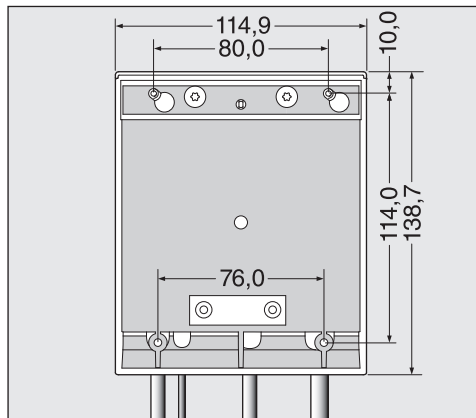**Wall mounted:**

- Using the drawing as a guide, fit the screws the correct distance apart (80.0 mm) and suspend the dispensing module from these.

**Surface mounted:**

Two slots 76.0 mm apart are provided at the bottom of the dispensing module for fixing the module to a thin carrier or plate. Screws 3.9 mm in diameter (not supplied) are required to secure the module.

- Drill two holes through the plate and screw the module firmly in place from underneath.

## en - Installation

---

**Installation inside an adjacent unit:** If more than one dispensing module is being used, or if the liquid agent canisters are to be stored with the module, a suitable unit is available from Miele.

Make sure hoses are laid without kinks.


## Connecting the dispensing modules

Please note that all dispensing modules and accessories supplied with them must only be fitted and commissioned by the Miele Customer Service Department or authorised service technicians.

The DOS G 80 module can be connected as standard to the washer-disinfectors and commercial dishwashers listed under “Areas of application for dispensing modules”.

The following accessories, which are supplied with the relevant dispensing module, must be used for connection to commercial dishwashers marked with an asterisk \*.

### DOS G 80 with hose adapter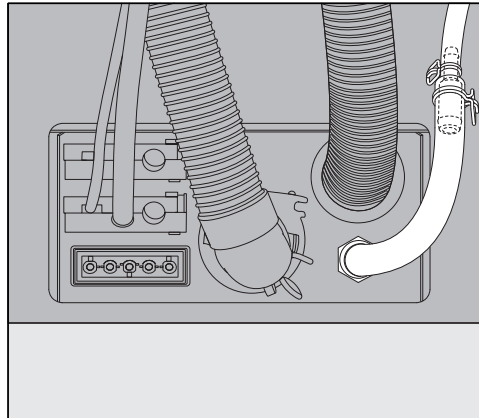


The hose adapter for dispenser connections with larger diameters is supplied with the DOS G 80 and must be used with G 8066 and PG 8067 commercial dishwashers.

Using the adapter increases the diameter of the dispensing hose to match the diameter of the connection.

### DOS G 80 ProfiLine with hose adapter and dispenser connection

The dispenser connection and the hose adapter are supplied with the DOS G 80 ProfiLine and must be used with ProfiLine commercial dishwashers.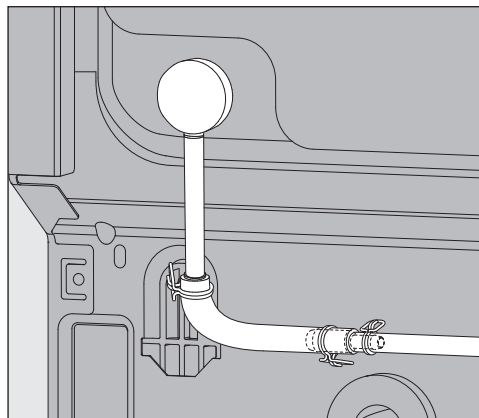

- Remove the blind stopper and fit the dispenser connection.

The dispenser connection may be hidden behind the back panel of the commercial dishwasher.

## en - Connection


- Unscrew the back panel and disengage the area around the dispenser connection.

### **When setting the dosage please note:**

For cleaning agent dosage the dispensing concentration must be doubled in the controls for PG 8080, PG 8081i, PG 8082 SCi and PG 8083 SCVi machines.

See the specifications for setting the dispensing concentration in the operating instructions.

### **Strain relief for dispensing hoses**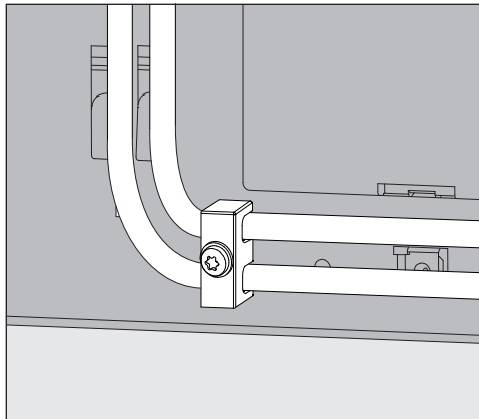


The strain relief provided with the dispensing modules for the dispensing hoses is fitted to the bottom right-hand side of the back of the machine to prevent the hoses from dislodging themselves from the dispenser connections during use.

### **Connecting the dispensing hoses**

The dispensing module is tested at the factory for functionality. Any residues of the harmless test medium are a result of this testing and not evidence of previous use.

- Pull the safety cap off the end of the hose and remove the protective foil from the siphon.
- Connect the hose with the dispenser connection to the cleaning machine.
- Connect the module to the machine's power supply.

Connections for the dispensing hoses are located on the back of the cleaning machine. These are described in detail in the operating instructions for your cleaning machine.

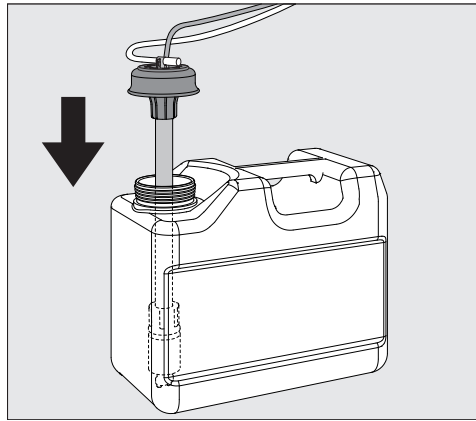- Fill or replace the container, fit the siphon and secure the lid.
- Feed the siphon into the canister until it reaches the bottom.

Follow the instructions for venting the dispensing system and setting the dosage in the operating instructions.

**Colour coding on the dispensing hoses**

To differentiate between the canisters when using more than one dispensing module, the hoses must be colour-coded using the colour labels supplied.

- Stick a label near the siphon and one of the same colour close to the free end of the hose to enable clear identification.

**Connecting the power supply**

The power supply connections are located on the back of the machine.

(They are described in detail in the operating instructions for your cleaning machine.)

- Connect the module to the machine's power supply.

**DOS G 80/72**

DOS G 80/72 is supplied without a plug. It is designed for direct connection to the PG 8172 tank dishwasher.

**Maintenance**

The dispensing system should be serviced at the same time as the cleaning machine by Miele Service.

## en - Technical data

Height	138.7 mm	
Width	114.9 mm	
Depth	114.2 mm	
Weight (net)	1.5 kg (incl. siphon, hose and cable)	
Connection	Connection to cleaning machine	
Voltage, Connected load, Fuse rating	200 - 230 V 50 Hz 12 W 16 A	200 - 240 V 60 Hz 10 W 16 A
Throughput (nominal)	55 ml/min	
Ambient temperature	5 °C to 40 °C	
Relative humidity maximum linear decreasing to	80 % for temperatures up to 31 °C 50 % for temperatures up to 40 °C	
Installation height	as per cleaning machine	
Pollution degree	2	
Safety class (according to DIN EN 60529)	IP21	
Test certificates awarded	EMC, VDE, radio and television suppressed	
CE-mark	2014/35/EC Low voltage directive	
Manufacturer's address	Miele & Cie. KG, Carl-Miele-Straße 29, 33332 Gütersloh, Germany	

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