
# Mobile injector unit E 350


## Items supplied:

1 water feed pipe	15 x E 351 jets, 4.0 x 160 mm
1 seal	15 x E 353 clips for 4.0 x 160 mm jets, consisting of holder + clamp
1 mobile unit E 350	18 x E 352 jets, 6.0 x 220 mm
1 jet for flushing out powder dispenser	18 x E 354 clips for 6.0 x 220 mm jets, consisting of holder + clamp

## Installation: Tools required: Spanner SW 9, cross-tip screwdriver

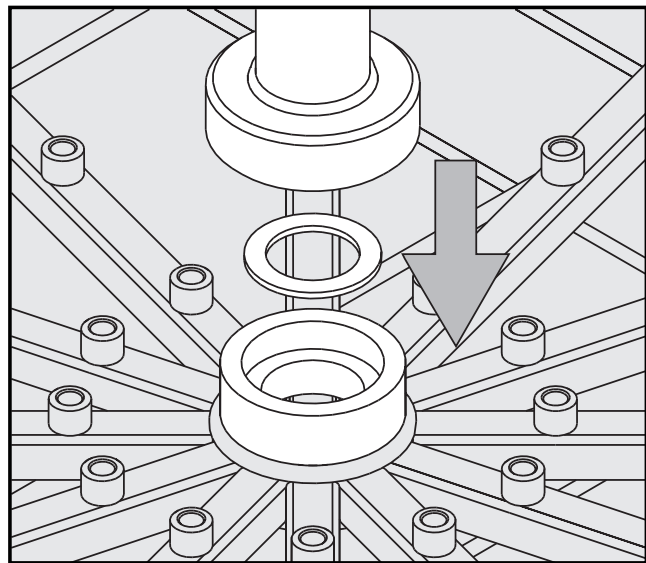
- Fit the seal in the holder in the mobile unit.
- Then fit the water feed pipe into the holder.
- Tighten the nut by hand.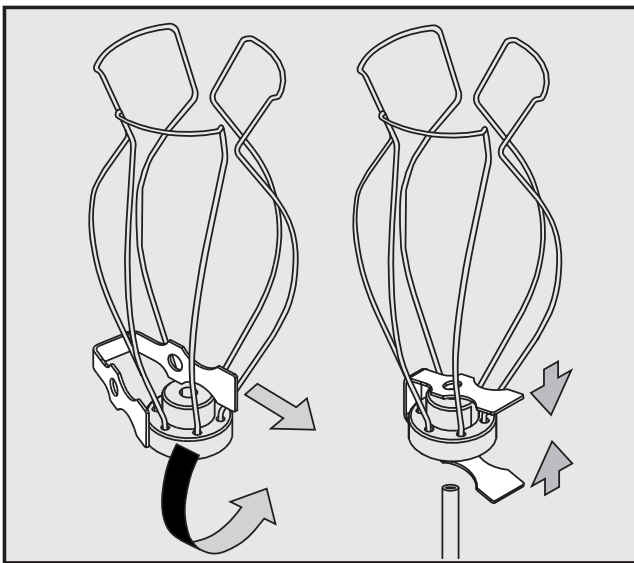


The mobile unit is supplied with two sizes of clamps, holders and jets. Make sure they are put together in the correct order.

- Push the clamp into the holder as shown above.
- Then squeeze the clamp together and feed the injector jet through the holes in the clamp.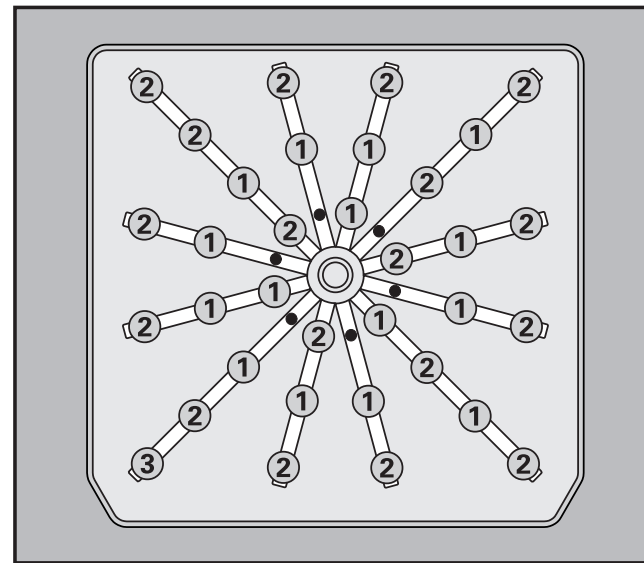

- Use the spanner to secure the jets into the positions required. The overview above gives an example of possible positions. Start from the centre and work your way outwards in ever increasing circles.
  - ① E 351 jets, 4.0 x 160 mm
  - ② E 352 jets, 6.0 x 220 mm
- Screw the jet for flushing out the powder dispenser in position ③ and secure it to the spray arm with a cross-tip screwdriver.

## Areas of application:

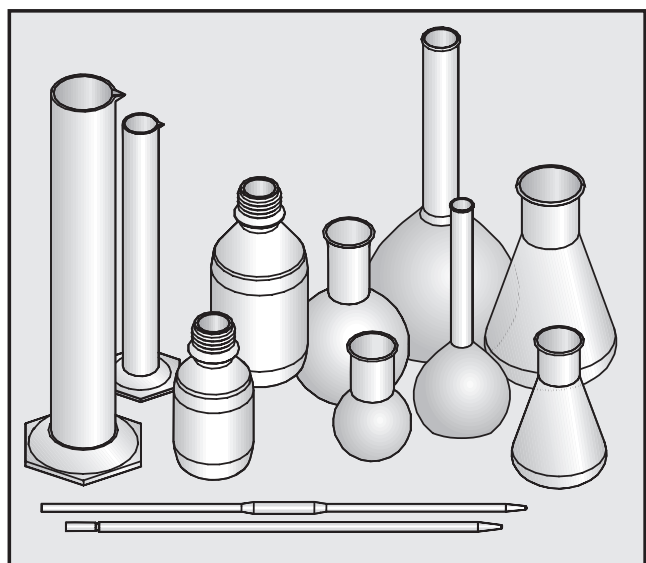
To ensure adequate water pressure all positions need to be equipped with jets.


- The mobile injector unit can hold up to 33 pieces of laboratory glassware.
- Different volume glassware can be processed, however, the maximum volume per piece is 1 litre.
- To process pipettes special sleeves are needed, MIBO E 336. They are available to order.
- Laboratory glassware, which because of its design cannot be flushed through properly, will not be sufficiently cleaned and disinfected in this mobile unit.
- This mobile unit should only be used for processing laboratory glassware which has been declared as suitable for machine processing by the glassware manufacturer.

The jet for flushing out the powder dispenser is designed to rinse out the powder dispenser in the machine door after use (if applicable to the machine the unit is used in). This jet must therefore not be used for cleaning glassware.


## Place the laboratory glassware over the jets:

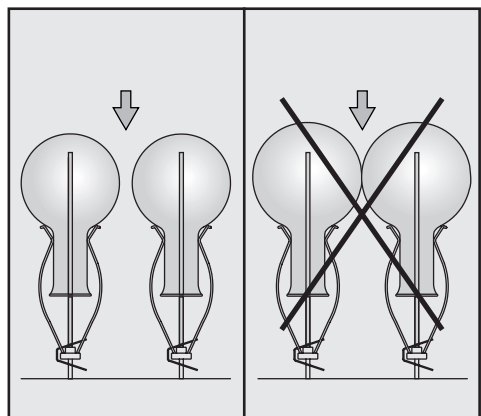
Do not let the individual pieces of glassware touch each other.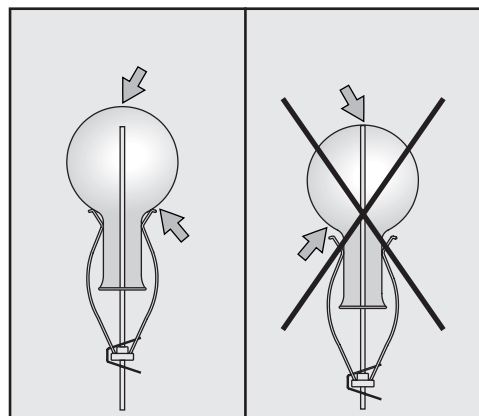

Push the clips upwards, if necessary, to prevent the tip of the jet from contacting the bottom of the glassware.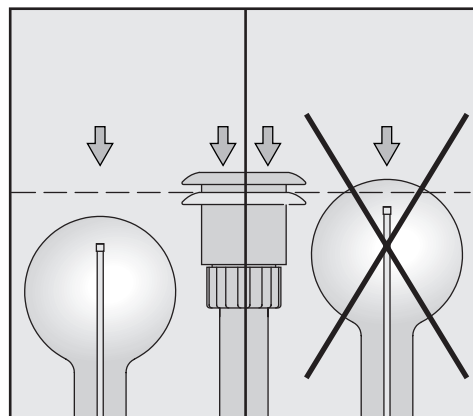

Glassware must not sit higher than the top of the water feed pipe (with the water feed adapter pressed downwards).

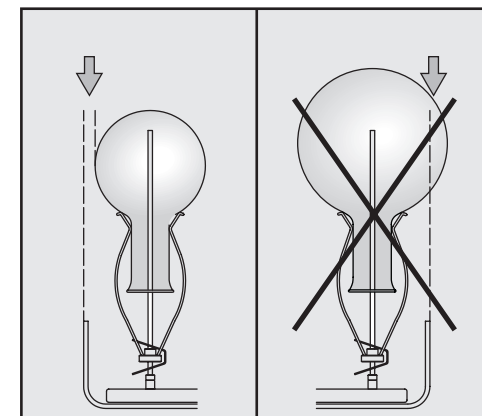Glassware must not extend beyond the sides of the mobile unit.

⚠ Carefully read the Warning and Safety instructions in the Operating Instructions for your washer-disinfector as well as any instructions provided by the glassware manufacturer.  
Any broken glassware can cause serious injury.  
All glassware should be properly inspected and tested for cleanliness. Do not rely on a visual check alone.

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