
Lower basket U 175/1

Items supplied: This lower basket is supplied complete. The following individual parts are included:

- | | |
|-------------------|--|
| 1 water feed pipe | 1 jet for flushing out powder dispenser |
| 1 seal | 33 x E 351 jets, 4.0 x 160 mm |
| 1 lower basket | 33 x E 353 clips for 4.0 x 160 mm jets, consisting of holder + clamp |


Installation: **Tools required:** Spanner SW 9, cross-tip screwdriver

The lower basket should be checked from time to time. Should any of the individual components have become loose or damaged in any way, they can be re-secured or replaced by following the instructions below.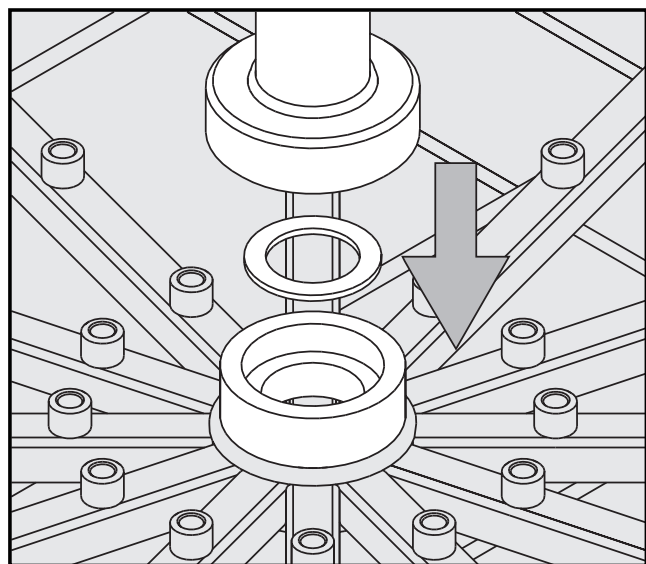


The water feed pipe will be easier to remove and refit if injector jets are first removed from the centre of the basket.

- Fit the seal in the holder in the basket.
- Then fit the water feed pipe into the holder.
- Tighten the nut by hand.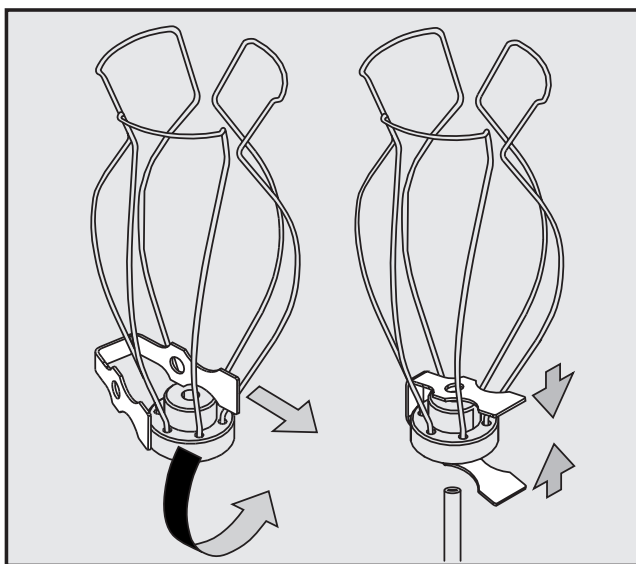


- Push the clamp into the holder as shown above.
- Then squeeze the clamp together and feed the injector jet through the holes in the clamp.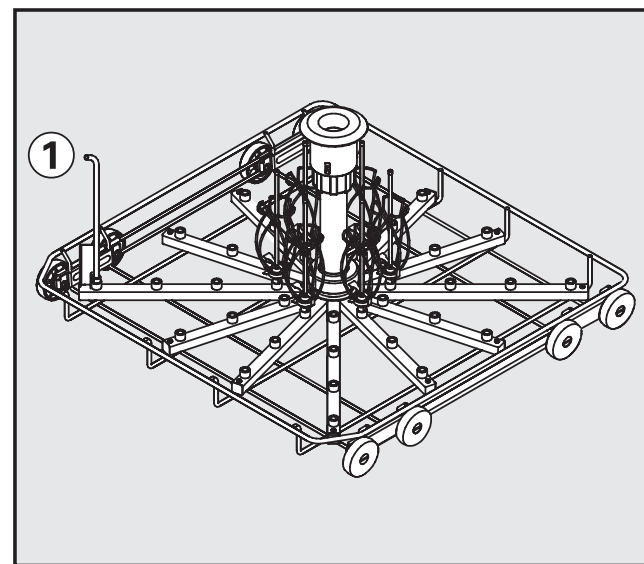

- Screw the injector jets into position using the spanner. Start from the centre and work your way outwards in ever increasing circles.
- Screw the jet for flushing out the powder dispenser in position ① and secure it to the spray arm with a cross-tip screwdriver.

To ensure adequate water pressure all positions need to be equipped with jets.


Areas of application: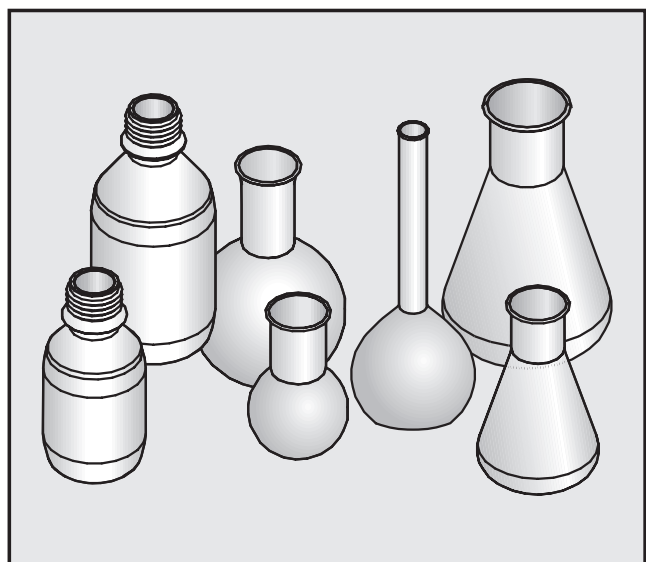


- Lower basket U 175/1 can only be used in conjunction with upper basket O 175 or O 187.
- This basket can be fitted with up to 33 pieces of laboratory glassware with a maximum volume of 200 ml or maximum height of 200 mm.
- Laboratory glassware, which because of its design cannot be flushed through properly, will not be sufficiently cleaned and disinfected in this mobile unit.
- This mobile unit should only be used for processing laboratory glassware which has been declared as suitable for machine processing by the glassware manufacturer.

The jet for flushing out the powder dispenser is designed to rinse out the powder dispenser in the machine door after use (if applicable to the machine the unit is used in). This jet must therefore not be used for cleaning glassware.


Place the laboratory glassware over the jets: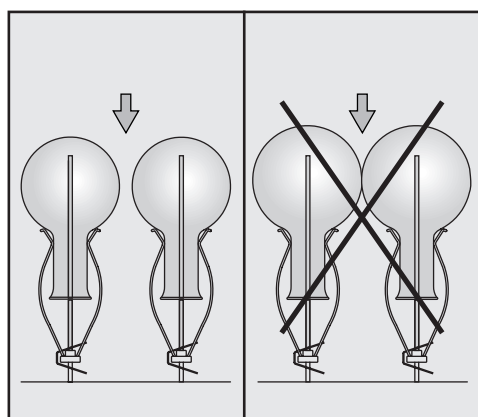


Do not let the individual pieces of glassware touch each other.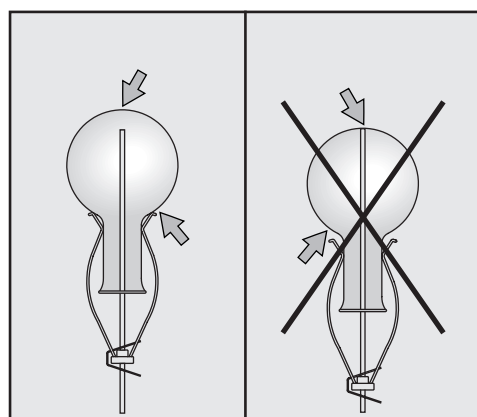

Push the clips upwards, if necessary, to prevent the tip of the jet from contacting the bottom of the glassware.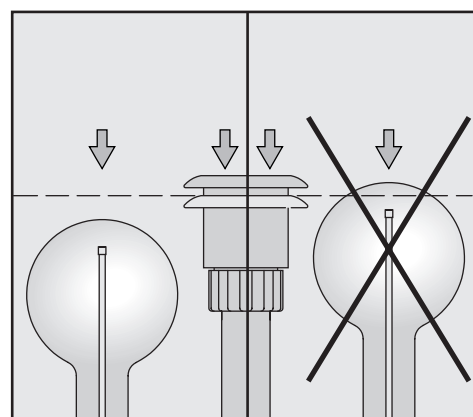

Glassware must not sit higher than the top of the water feed pipe (with the water feed adapter pressed downwards).

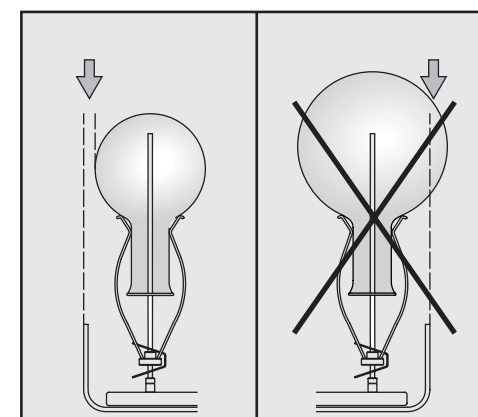Glassware must not extend beyond the sides of the mobile basket.

⚠ Carefully read the Warning and Safety instructions in the Operating Instructions for your washer-disinfector as well as any instructions provided by the glassware manufacturer.

Any broken glassware can cause serious injury.

All glassware should be properly inspected and tested for cleanliness. Do not rely on a visual check alone.

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